

InSpec™ Biocides

Technical Data of InSpec Products by Chemistry

Aqueous High Level Disinfectant



InSpec™ AN

Sterile Diamine, Disinfectant Cleaner

Low Residue, Quat Free Disinfectant/Cleaner, Broad Spectrum Disinfectant

Use Instructions

InSpec™ AN is designed for spraying, wiping and mopping applications. Hold spray approximately 1.5cm to 20cm from area to be treated. 5L versions, pour into an appropriate container for mopping. 100ml concentrate added to 2.5L of water is required for mopping. Apply to surface to ensure complete coverage.

Technical Data

Alcohol, chlorine and QUAT-free disinfectant cleaner. 4% RTU Solution, pH 8.8-9.2, Concentrate pH 10.0.

Storage & Safe Handling

Store upright in original closed containers, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Safety Data Sheet (SDS).

Efficacy

See summary table in Section 1 of this brochure.

Material Compatibility

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical Dossier.

Benefits

- An exceptional non-oxidising biocide
- Bactericidal and Fungicidal
- Clean and disinfects
- 3 months validated 'in-use' shelf life
- Contents protected using 'bag-in-bottle' presentation with premium range
- 1L and 5L formats manufactured with WFI quality water
- Multiple bags for cleanroom transfer
- Ideal rotational partner for InSpec QT, HA and OX
- Manufactured in accordance with GMP

Technical User Information

InSpec™ AN is a sterile disinfectant in pre-diluted and concentrated formats. InSpec™ AN is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.

The solution is filtered through a 0.2 micron filter and gamma irradiated to a validated process (at 25-45kGy) to give a sterility assurance level SAL of 10⁻⁶.

Trigger sprays supplied with 900ml solution in a 1L container.

InSpec™ AN is supplied with Certificates of Conformity, Analysis, Irradiation and Sterility. All formats of AN have a stabilised 24 month shelf life.

For screw cap bottles use the entire contents after opening.

InSpec™ AN Sterile Products		
Product	Sterile Part No.	No. per case
InSpec AN 900MLS PREMIUM	ANWFI20-1LS	6
InSpec AN 5L	ANWFI20-5LS	2
InSpec AN Concentrate	ANCNC20-100	50
InSpec AN Burstable Pouch - 15 wipes	ANWFI20BP15S	20

InSpec™ AN Non-Sterile Products	
Non-Sterile Part No.	No. per case
NSANWFI-1L	6
NSANWFI-5L	2
NSANCONC-100ML	50
N/A	N/A

Wipe Media Type	Individual Wipe size (mm)
N/A	N/A
N/A	N/A
N/A	N/A
Polycellulose 68gsm	230 x 230