

**InSpec™ Biocides**

# Technical Data of InSpec Products by Chemistry

## Aqueous Sporicidal Agents



# InSpec™ OX

## Sterile Sporicidal Preparation Hydrogen Peroxide 6% v/v and Peracetic Acid 0.4% v/v

Rapid action - two minute kill time. Achieves the EN sporicidal standard - 3 log reduction in 2 minutes

**Use Instructions**

InSpec™ OX is designed for spraying, wiping and mopping applications. Hold spray approximately 15cm to 20cm from area to be treated. 5L versions, pour into an appropriate container for mopping.

**Technical Data**

Appearance: Colourless clear liquid with a characteristic strong acetic odour. pH 1.4.

**Storage & Safe Handling**

Store upright in original closed containers, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Safety Data Sheet (SDS).

**Efficacy**

See summary table in Section 1 of this brochure.

**Material Compatibility**

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical Dossier.

**Benefits**

- One of the most effective sporicides available on the market
- All formats classed as an irritant.
- Low Hazard
- Stabilised solution - 24 months shelf life
- 3 months validated 'in-use' shelf life
- Contents protected using 'bag-in-bottle' presentation with premium range
- 1L and 5L formats available as Ready to Use
- Multiple bags for cleanroom transfer
- Ideal rotational partner for InSpec QT and AN
- Manufactured in accordance with GMP

**Technical User Information**

InSpec™ OX is a sterile sporicide available ready to use.

InSpec™ OX is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.

The solution is filtered through a 0.2 micron filter and aseptically filled into pre-gamma irradiated packaging (25-45kGy) and tested to give a sterility assurance level SAL of 10<sup>-6</sup>.

InSpec™ OX is supplied with Certificates of Conformity, Analysis, Irradiation and Sterility. All formats of OX have a stabilised 24 month shelf life.

For screw cap bottles use the entire contents after opening.

InSpec™ OX Sterile Products		
Product	Sterile Part No.	No. per case
InSpec OX 1L PREMIUM Trigger Spray	OXPW40-1LS	6
InSpec OX 5L	OXPW40-5LS	2
InSpec OX Burstable Pouch - 15 wipes	OXPW40BP15S	20
InSpec OX Burstable Mop - 2 EF40 Mops	Z116004	5

InSpec™ OX Non-Sterile Products	
Non-Sterile Part No.	No. per case
NSIOXPW-1L	6
NSOXPW-5L	2
N/A	N/A
Z116003	5

Wipe Media Type	Individual Wipe size (mm)
N/A	N/A
N/A	N/A
N/A	N/A
100% Polyester microfibre	400 x 100